

Calibrating the Con-6

AUTOMATIC CALIBRATION:

- * Set meter to OFF.
- * Pour just enough 12,880 calibrating solution into test tube to cover the probe.
- * Insert probe into 12,880 calibrating solution.
- * Press [ON].
- * Press [CAL].
- * Press and hold [HOLD/ENTER] until reading is 12.88 or close.
- * Remove probe and rinse with distilled water.

MANUAL CALIBRATION:

- * Set meter to OFF.
- * Pour just enough 12,880 calibrating solution into test tube to cover the probe.
- * Insert probe into 12,880 calibrating solution.
- * Press [ON].
- * Press [CAL].
- * Use UP/DOWN keys to adjust display to calibrating solution.
- * Press [HOLD/ENTER] to confirm.
- * Remove probe and rinse with distilled water.